
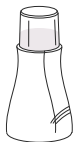

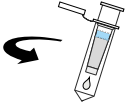
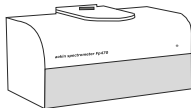


Sample preparation with **aokin** QuickClean

Extraction	1. Weighing out		directly into extraction beaker 5 g sample + 32 g aokin mycontrol DON Extraction solvent (31.8 ml at 20 °C)	~ 1 min
	2. Extraction		mix 1 minute	~ 1 min
	3. Filtration		collect filtrate (discard filter cake)	~ 1 min
Purification	4. Quick clean		400 µl filtrate to aokin QuickClean column centrifuge 2 minutes at 5000 x g	~ 2 min
Measurement	5. Measurement		aokin spectrometer FP470	~ 3 min